

# Genesis Quick Start Guide

## **Charger**

Pull and insert into a wall plug (110VAC)

Before shipment, insert cord and wall adapter back into charger cavity.

## **On-Off Switch**

Lockable tamper-resistant switch

0 -off / I -on  
Rotate Clockwise to turn on.


## **Payload Canister**

Insert vials into openings.

Insert foam plug to secure vials in place.

## **Data Port**

Shipper programming & data retrieval port.

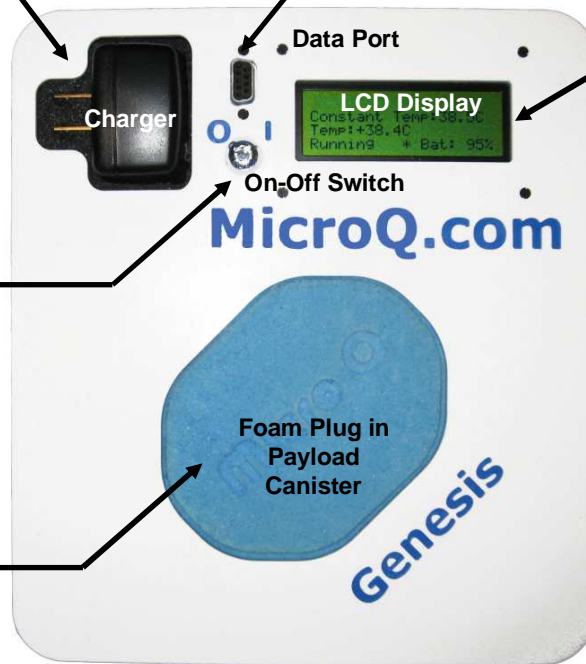

## **LCD Display**

Follow instructions on screen.  
Screen display may vary. See examples

## **Start up screen**

```
Micro Q LLC
Genesis
Version 1.1
Copyright 2009
```

## **Running screen**

```
User Name
33 to 180, 21.0 Hr
Time:58 T:+33.0 C
Unplugged Bat:100 %
```

## **Information Screen**

```
Low Battery Power
DO NOT SHIP
PLUG IN Charger
Wait for Full Charge
```

microQ.com

# Genesis Quick Start Guide

## **Instructions**

### **Shipment.**

1. Charge device fully.
2. Unplug device.
3. Turn on device to initiate temperature protocol.
4. Plug charger into outlet. Charge device until just prior to shipment.
5. Pre-condition canister to the appropriate temperature.
6. Load samples
7. Insert Genesis into exterior ventilated box
8. Just prior to shipment, unplug and stow charger and cord.
9. Tape box closed making sure to not block ventilation openings.
10. Attach shipping label to the top of box.

### **Receiving.**

1. Open exterior box and check battery condition.
2. Gently, remove charger and plug into outlet.
3. Grasp foam plug in center of faceplate and pull to reveal sample containment area.



microQ.com